

DRV-Z1 Near-Coaxial Illuminator - Instructions

The DRV Near-Coaxial Illuminator will work with DRV-Z1 zooms DRV200 and DRV201.

Fitting

- Remove any fitted DRV-Z1 objective lens.
- Fit the objective lens to the illuminator.
- Fit the illuminator to the objective mount of the DRV-Z1.
- Connect the power cable to the supply.

A video of the installation process can be accessed here:

<https://www.youtube.com/watch?v=-UoBFuv7X9E>

Usage

- Power on the illuminator
- Use the DRV-Z1 keypad to turn off the DRV-Z1 ring-light.
- Use the Near-Coaxial controller to adjust the brightness.